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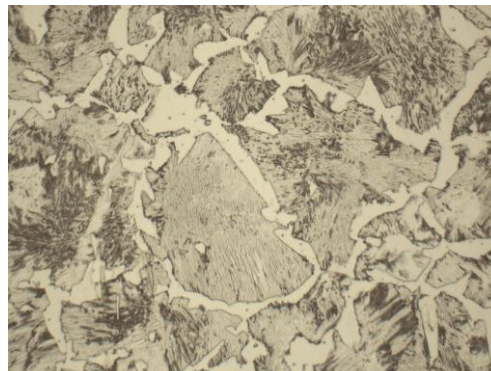
 **Struers**

Pederstrupvej 84, DK-2750 Ballerup

Transcopy Replica Foil creates a permanent replication of microstructures. The metallisation of the back side makes **Transcopy Replica Foils** highly reflective and gives a high contrast of strongly etched surface detail.

NEW colourless type!

Instruction for use

- The surface to be replicated must be polished and etched. It is recommended to over etch the area of interest.
- *Using the pipette:*
 - Apply 2 drops of Transcopy Liquid onto the Transcopy Replica Foil.
 - Let the liquid flow around on the Transcopy Replica Foil so it is evenly distributed (avoid wetting the self-adhesive backing).
- *Using the spray nozzle:*
 - Hold the bottle approx. 15 cm / 6" from the area to be wetted and spray twice, directly onto the sample surface or on the Transcopy Replica Foil.
- Wait a few seconds.
- Press the foil hard against the etched area with a thumb for 30 seconds.
- Wait 4-5 minutes then carefully remove the Transcopy Replica Foil.
- Remove the backing paper and adhere the Transcopy Replica Foil to a microscope slide.

Tips

- The contrast can be increased on larger optical microscopes using interferential interference contrast (DIC) at high magnifications or filters (e.g. grey or green ones).
- Replicas can be stored for decades and subsequently used for comparative analyses, provided they are protected against dust, heat and humidity.

Transcopy Liquid Safety Data

Please download a copy of the **Safety Data Sheet** from the Library section on www.Struers.com and read carefully before use. Access the Struers Library from the top menu on the website.

Please check the Library periodically for updates to the Safety Data information.

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